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INSTRUCTION MANUAL

STEREO ZOOM
MICROSCOPE XTS2021


Thank you for purchasing the Stereo Zoom Microscope.
Please read this instruction manual thoroughly before use to insure proper operation and long service.

■NOTES

- This is a precision instrument – Handle With Care. Use caution to avoid shock or excessive forces on instrument when in use or transport.
- Avoid direct sunlight or areas with high temperatures or high levels of dust. Keep in location where temperate and humidity are stable.
- Keep lens surface clean from dirt and fingerprints.
- Clean the lens gently with lens cloth or soft cotton cloth.
- Do not turn the left and right Zoom Controls in opposite directions; this will damage the instrument.
- When not in use store with protective cover, in location that has low moisture in order to prevent internal mold growth.
- Adjust the Light Power Dial to adjust the brightness of the epi-illuminator for reflected light.
When first turned on, rotate the dial to check that the power to the light is on.
- Lighting unit not certified for use outside of Japan.

■SPECIFICATIONS

Rated Voltage	AC100V 50/60Hz
Magnification	6.5~45× (for 10× eyepiece) 13~90× (for 20× eyepiece)
Objective Lens	0.65~4.5× (Zoom Ratio 1:7)
Eye-piece Lens	WF10× (FOV = 20mm)・WF20× (FOV = 10mm)
Working Distance	95mm
Eyepiece Angle	45°
Focus Adjustment	Movable Body
Stand	Rack and Pinion Type
Stage Plate	φ95mm×2pcs ①Ground Glass ②black/white (plastic)
Eye Width Adj. Range	55~75mm
Diopter Adj. Range	±5diopter
Illumination	Epi-illuminator: 6V・15W Halogen bulb Transmitted: 5W ring light
Base	L245×W190×H60mm
Weight	approx.5.7kg

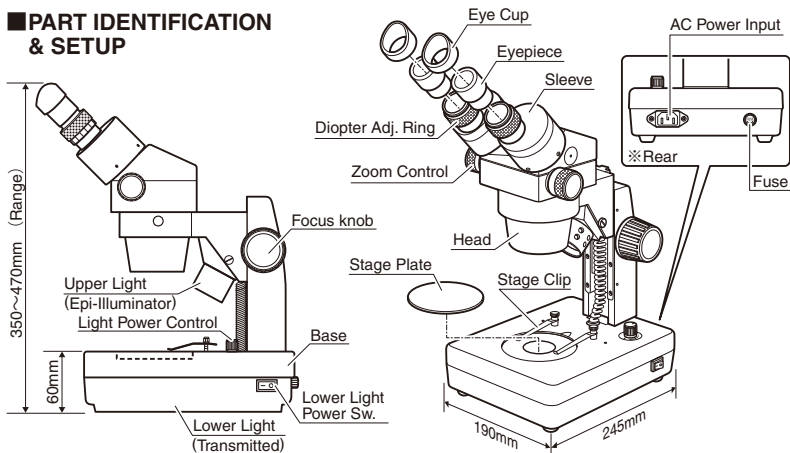■PART IDENTIFICATION
& SETUP

■OPERATION

SAMPLE
PLACEMENT

- ① Place the sample in the center of the Stage Plate
- ② If required, use Stage Clips to hold sample in place.

EYE CUP
USE

- For use without eyeglasses : Hold Eyepiece and rotate the Diopter Adjustment Ring to adjust diopter setting. Align the Eye cups to point toward user and look through the Eyepiece with eyes brought close.
- For use with eyeglasses : Remove Eye cups to use.

EYE WIDTH
ADJ.

- Hold both sleeves in each hand and look through the Eyepieces. Adjust the Sleeves to align the left and right field of view seen through the Eyepieces.

FOCUS &
DIOPTRER
ADJUSTMENT

- ① Rotate the Zoom Control for highest magnification.
- ② Look through the right eyepiece with the right eye and adjust the focus using the Focus Knob.
- ③ Rotate the Zoom Knob for lowest magnification.
- ④ Using the Right Diopter adjustment, adjust the focus if required.
- ⑤ Now rotate the Zoom Control back to highest magnification and check the focus.
Repeat steps ② ~ ④ to achieve the best focus accuracy.
- ⑥ Rotate the Zoom Control for highest magnification. Look through the left Eyepiece with the left eye, and adjust focus using the Left Diopter Adjustment Ring.

FUSE
REPLACEMENT

- In case of fuse failure, it can be replaced with a commercially available 250V 2A fuse.

SC Niigata seiki Co., Ltd.
5-3-14, Tsukanome, Sanjo, Niigata, Japan, 955-0055
Tel. : +81-256-33-5522 Fax. : +81-256-33-5518
MAIL intl.sales@niigataseiki.co.jp
URL <http://www.niigataseiki.co.jp>